

SUMMIT MEDICAL

Equipment Company



SLIDING TOP RODENT INDUCTION CHAMBERS

OPERATING INSTRUCTIONS

PN: AS-01-0530-LG RAT & MOUSE

PN: AS-01-0530-SM MOUSE

DIMENSIONS: 5.25" H X 5.375" W X 11.25" F/B

4.625" H X 4.25 W X 9.25" F/B

VOLUME: 5.4 liters

3.0 liters

1. Place the chamber on a flat surface with the opening away from the operator, so that the sliding top slides toward the operator.
2. When the animal is fully anesthetized turn the Isoflurane off, and allow about five seconds for Oxygen flow to help reduce Isoflurane concentrations in the chamber.
3. Lift the chamber at the end closest to the operator and gently roll the animal down to the far side.
4. Placing your hand over the far end of the chamber slide the top back just far enough to insert your hand and lift the animal out. Then immediately close the top.
5. By keeping the opening as far away from the operator as is reasonably possible, and reducing the time that the chamber is open as much as possible, the risk of exposure to the operator will be reduced.